




CAVACam

An Ultra-Clean Camera for UCV Vacuum Applications

Introduction

CAVACam is a customisable CAmera for VAcuum applications, specifically designed to be compliant with ultra-clean vacuum (UCV) standards. The CAVACam has been manufactured using vacuum-compatible materials which are low-outgassing and minimise contamination risk. Available with a wide range of detectors, the modular design can be easily adapted to suit a range of scientific applications.

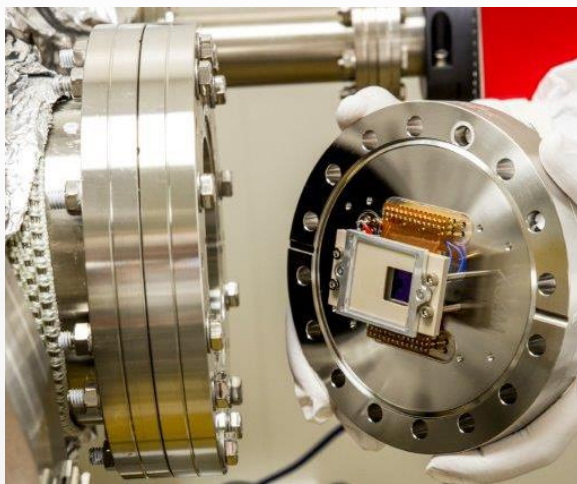

CAVACam (F-model)

The CAVACam is available as a conflat flange model (F-model) or as a hermetically sealed, fully vacuum-immersible model (S-model). Detectors are available in both front-illuminated (FI) and back-illuminated (BI) formats to cater for all imaging requirements. BI detectors may also be available with no AR coating for improved soft X-ray performance. EMCCDs are available for low-light imaging and single photon detection applications.

XCAM has a wealth of experience and expertise manufacturing cameras for ultra clean vacuum applications having successfully developed vacuum cameras for many prestigious scientific and industrial institutions around the world.

Key Features


- Clean construction to UHV, XHV or UCV standards as required
- Guaranteed not to contaminate your vacuum chamber if operating instructions are properly followed
- Available with a wide range of EMCCD and CCD detectors, front- or back-illuminated
- Electronic drive systems suitable for:
 - high flux and high frame rate with exceptional linearity
 - extremely low noise levels at slow speeds for sensor-limited noise performance
 - single photon detection and sub-pixel spatial resolution, as demonstrated by the RIXSCam, one of XCAM's signature advanced camera systems
- Multiple camera heads and/or detectors can be operated synchronously
- Systems are individually optimised for each detector type to give the best performance
- Water-cooled Peltier capable of cooling down to -50°C if required
- Cryocooled solutions also available capable of cooling to -110°C
- Modular design enables easy customisation to cater for individual requirements
- RGA scan can also be conducted if required



CAVACam

An Ultra-Clean Camera for UCV Vacuum Applications


CAVACam F-Model (Conflat Flange)

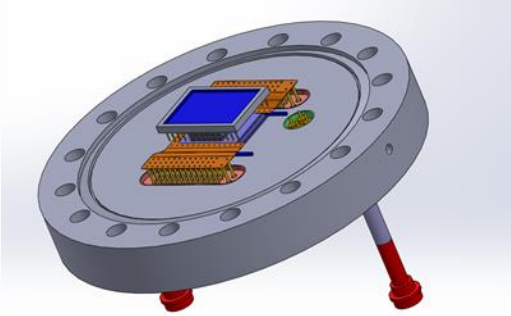

The CAVACam F-model is a camera head which is built on a conflat-style flange. The detector and cooling system are on the vacuum-side of the flange, with the electronics on the air-side of the flange. Hermetic electrical feedthroughs maintain the vacuum chamber integrity and ensure that the non-UCV compatible electronics remain safely away from your vacuum experiment.

CAVACam S-Model (Fully Vacuum-Immersible)

The CAVACam S-model is a camera head which can be entirely immersed inside an experimental chamber. Hermetically sealed in a stainless steel 'box', the non UCV-compatible electronics are safely sealed away inside the vacuum to prevent contamination. The detector is on the outside of the box and fully immersed in the vacuum chamber. The S-model is ideal for applications where the detector needs to be fully immersed inside the vacuum chamber and an F-model CaVACam may not be possible due to spatial limitations of the experimental set-up.

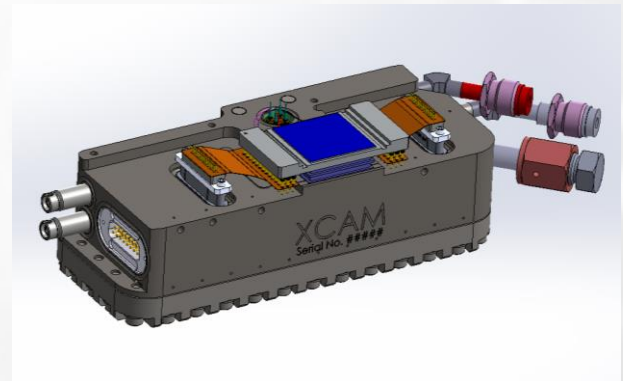

CAVACam Detector Range

The 2008, 2010 and 2011 series have been designed to cater for a variety of imaging requirements. Each CAVACam series is available in both the F-model and S-model. A wide range of detectors, in both full frame and frame transfer, are available within each series; XCAM have special expertise in selecting the ideal detector for the job.

- | | |
|--------------------|--|
| 2008 Series | <ul style="list-style-type: none">• <i>Extremely low noise at slower readout speeds</i>• Up to 4 channels/output nodes for readout speeds of up to 180KHz (lowest possible readout noise)• Up to 2 channels/output nodes for slightly higher readout speeds of 1-2 MHz |
| 2010 Series | <ul style="list-style-type: none">• <i>Ideal for high flux, high frame rate applications</i>• 4-channel operation at up to 3 MHz per output node (5 MHz coming soon)• Excellent linearity over whole range |
| 2011 Series | <ul style="list-style-type: none">• <i>Ideal for low-light applications where single photon detection may be required</i>• Individual photon-centroiding to within a fifth of a pixel for relevant applications• 4-channel operation at up to 3 MHz per output node (5 MHz coming soon)• Excellent linearity over whole range |

© 2017 XCAM. No part of this publication may be reproduced without prior permission in writing from XCAM. Whilst XCAM will endeavour to ensure that any data contained in this product information is correct, XCAM do not warrant its accuracy or accept liability for any reliance on it. XCAM reserve the right to change the specification of the products and descriptions in this data sheet without notice. Prior to ordering products please check with XCAM for current specification details. This product may be protected by patent. All brands and product names are acknowledged and may be trademarks or registered trademarks of their respective holders.

XCAM Ltd.
2 Stone Circle Road
Northampton
NN3 8RF
UK

Tel: 44 (0)1604 673700
Fax: 44 (0)1604 671584
Web: www.xcam.co.uk
Email: sales@xcam.co.uk